



Sensor Plug Assembly



CAUTION


The Pt100 of the sensor plug is made of glass and can break.
Do not bend or drop.

Follow these instructions to assemble the sensor plug, otherwise the glass probe may break.



1


Carefully unpack all parts of the sensor plug.

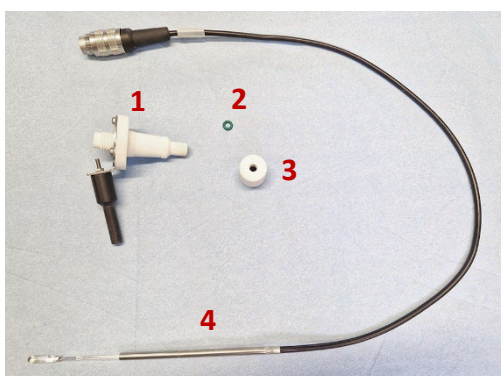

- 1 Distillation flask plug
- 2 O-ring
- 3 Union nut
- 4 Temperature sensor with glass Pt100

2


Slide the union nut **1** over the glass Pt100.
Then slide the O-ring **2** over the glass Pt100.

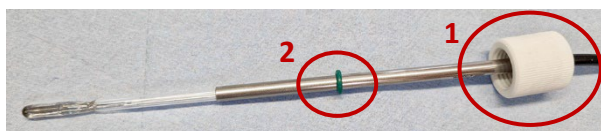

3

Carefully insert the glass Pt100 into the distillation flask plug.

When the Pt100 is in place, tighten the union nut.

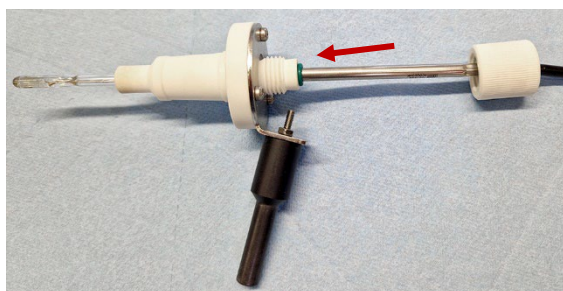

To adjust the height of the Pt100 in the distillation flask, loosen the union nut **1** and carefully pull or push the Pt100 to the desired position, then tighten the union nut again.

