

1 Heater Unit Replacement

If necessary, the replacement of the heater unit can be carried out directly by the user.

Required tools:

- Long screwdriver (optional)
- Long small pliers (optional)

1.1 Removing the Heater Unit

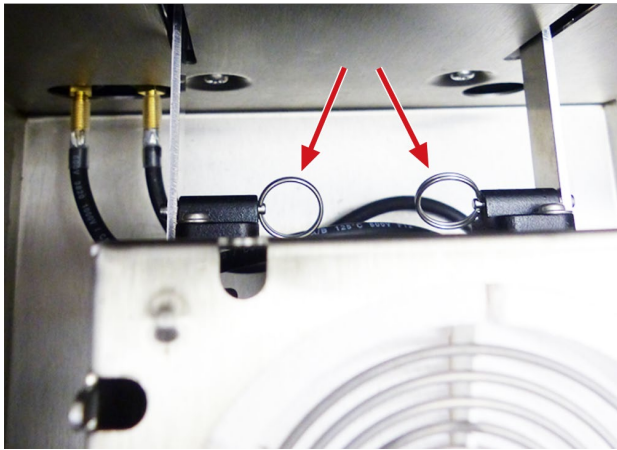


Fig. 1 Unlocking the heater unit

1. To remove the heater unit from the heater chamber you must pull two pins that are located just behind the heater unit. If you cannot reach them with your hand, you can use a screwdriver to pull them out.
2. The heater unit is also held in place by two hooks underneath it; to remove the heater unit, tilt it slightly towards you and push it back.

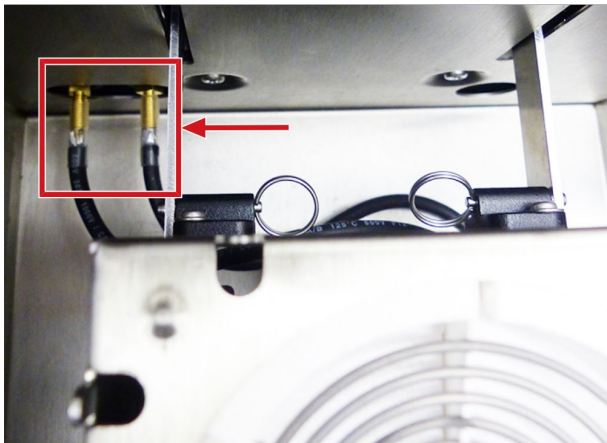


Fig. 2 Disconnecting the heater unit

3. Once the heater is detached, you can unplug the heater wires and remove the heater completely from the heater chamber.

1.2 Mounting the Heater Unit

To mount the new heater unit, follow the steps listed above in reverse order.

IMPORTANT:

Please check if the heater unit is locked by the two hooks underneath.

Contact us at

Anton Paar ProveTec GmbH

E-Mail: info.provetec@anton-paar.com

Website: www.anton-paar.com