

Setup Circulator Bath CD-B27 Set (mat. no. 180792, 180938, 188545)

Set up the connections of the circulator bath as shown in *Figure 1*.

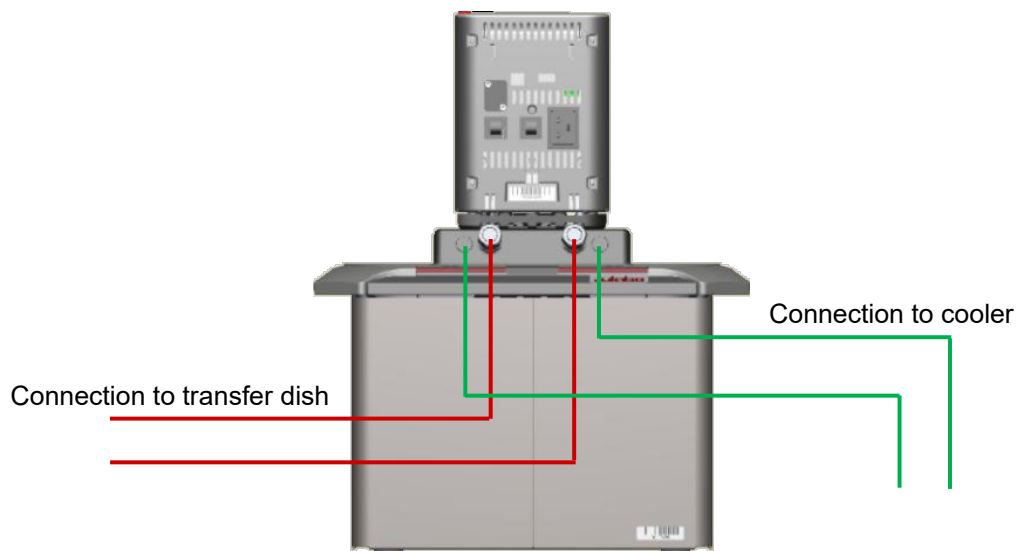


Figure 1: Back view of the circulator bath

Connect transfer dish, sample bath and cooler as shown in *Figure 2*.

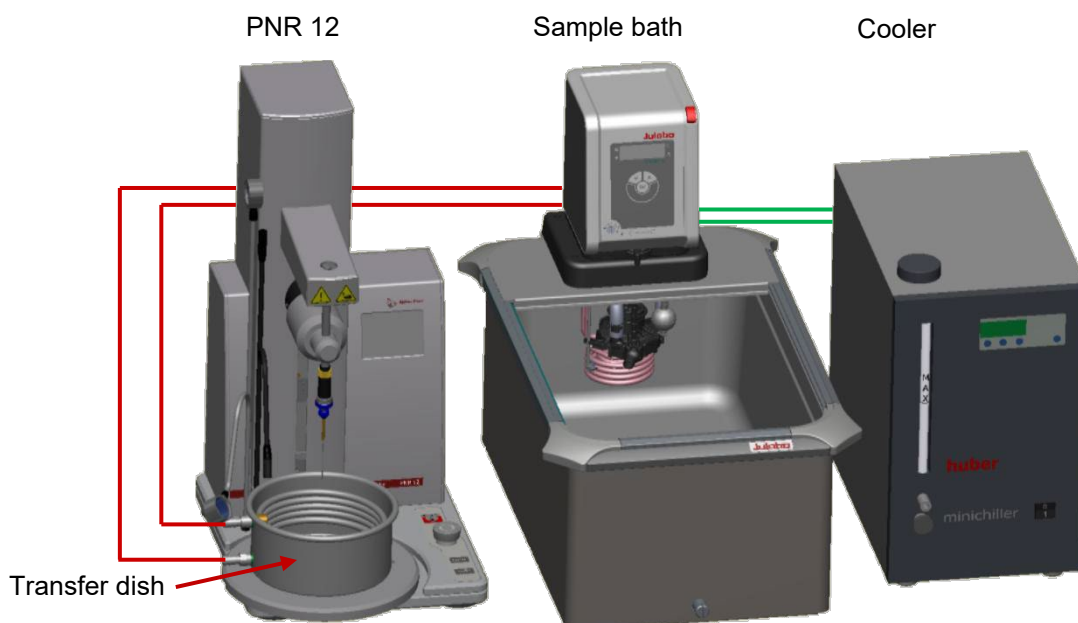


Figure 2: Setup of transfer dish, circulator bath and cooler

The cooler temperature should be set 5 °C lower than the sample bath temperature.

For example: bath temperature = 20 °C; cooler temperature = 15 °C

Contact us at

Anton Paar ProveTec GmbH

E-Mail: info.provetec@anton-paar.com

Website: www.anton-paar.com