


# A Chemist's Guide to Microwave Synthesis

Basics, Equipment & Application Examples

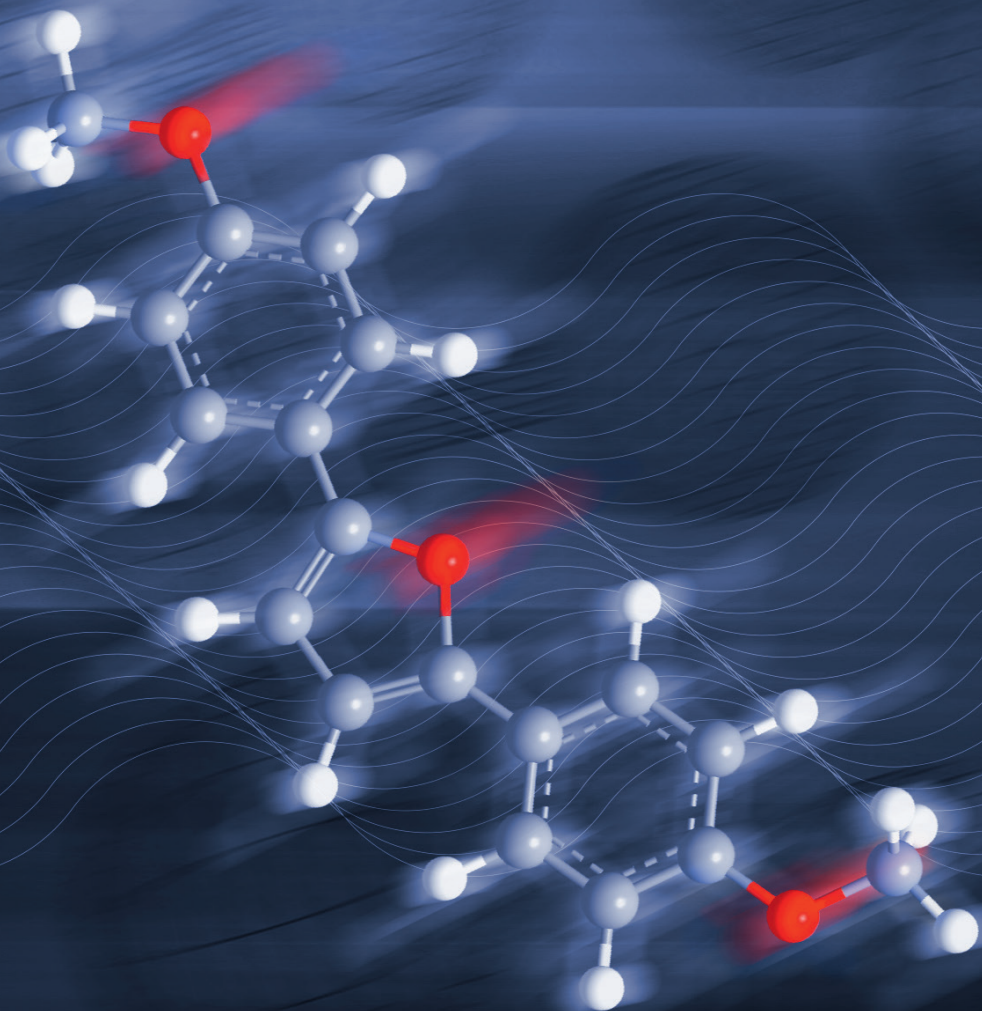

Jennifer M. Kremsner, Alexander Stadler

3<sup>rd</sup> Edition

# A Chemist's Guide to Microwave Synthesis

---

Basics,  
Equipment &  
Application Examples

by  
Dr. Jennifer M. Kremsner  
Dr. Alexander Stadler

- Copyright©2018 by Anton Paar GmbH, Austria.

All rights reserved. No part of this publication may be reproduced, stored in a retrieval system or transmitted in any form by any means electronic, mechanical, photocopying or otherwise without first obtaining written permission of the copyright owner.

- Published by **Anton Paar** GmbH. Printed in Austria.

**Anton Paar GmbH**

Anton-Paar-Str. 20

A-8054 Graz

Austria - Europe

Tel.: +43 316 257-0

Fax: +43 316 257-257

E-Mail: [info@anton-paar.com](mailto:info@anton-paar.com)

Web: [www.anton-paar.com](http://www.anton-paar.com)

*3<sup>rd</sup> Edition, January 2018*

# Contents

<b>1.</b>	<b>Introduction</b>	<b>6</b>
1.1.	About the Book	6
1.2.	General Aspects	6
1.3.	History of Microwave Synthesis	7
1.4.	Anton Paar and Microwave Synthesis	9
<b>2.</b>	<b>Principles of Microwave Heating</b>	<b>13</b>
2.1.	The Electromagnetic Spectrum	13
2.2.	Microwave Dielectric Heating	14
2.3.	Dielectric Properties	15
<b>3.</b>	<b>Benefits of Microwave Synthesis</b>	<b>18</b>
3.1.	Microwave vs. Conventional Heating	18
3.2.	Domestic Microwave Ovens vs. Dedicated Instrumentation	20
3.2.1.	<i>Safety</i>	20
3.2.2.	<i>Possibility of Superheating</i>	20
3.2.3.	<i>Excellent Parameter Control</i>	22
3.2.4.	<i>Automation and Parallel Synthesis</i>	23
3.2.5.	<i>Stirring</i>	23
3.2.6.	<i>Continuous Power Output</i>	24
3.2.7.	<i>Software Control</i>	25
3.3.	Controversial Statements and Features	27
3.3.1.	<i>Thermal and Non-thermal Microwave Effects</i>	27
3.3.2.	<i>Open Vessels vs. Closed Vessels</i>	28
3.3.3.	<i>Heating-while-cooling</i>	31
3.3.4.	<i>Power-controlled Reactions</i>	34

<b>4.</b>	<b>Equipment</b>	<b>36</b>
4.1.	Introduction	36
4.1.1.	<i>Multimode vs. Monomode</i>	36
4.1.2.	<i>Parallel vs. Sequential Processing</i>	38
4.1.3.	<i>The Importance of Accurate Temperature Measurement</i>	40
4.1.4.	<i>Silicon Carbide – Microwave Heating of Non-absorbing Reaction Mixtures</i>	43
4.2.	Monomode Reactors: Monowave Series	46
4.3.	Multimode Reactor: Masterwave BTR	49
4.4.	Multimode Reaction Platform: Multiwave PRO	51
4.4.1.	<i>High-throughput Accessory</i>	52
4.4.2.	<i>Parallel Synthesis and Parallel Scale-up</i>	54
4.4.3.	<i>Special Accessoires</i>	55
4.5.	Conventionally Heated Sealed-vessel Reactor: Monowave 50	56
4.5.1.	<i>Facts and Features</i>	56
4.5.2.	<i>Heating Efficiency</i>	58
4.5.3.	<i>Applications Examples</i>	59
4.6.	Summary – The Anton Paar Synthesis Family Philosophy	61
<b>5.</b>	<b>Getting Started – Frequently Asked Questions</b>	<b>63</b>
5.1.	How do I Find a Proper Starting Point?	63
5.2.	How do I Choose the Right Equipment?	65
5.3.	What are Typical Fields of Application?	68
5.4.	Is it possible to perform enantioselective reactions in microwave reactors?	69
5.5.	Which solvents can be used in microwave synthesis?	69
5.6.	How important is stirring in a microwave reactor?	70
5.7.	Can metals be safely heated in a microwave reactor?	70
5.8.	Can reactions which require inert gas atmosphere be performed in microwave reactors?	71

5.9.	Can solids be heated in a microwave reactor?	72
5.10.	Do I always need simultaneous internal and IR temperature measurement?	72
5.11.	Are there any known radiation hazards associated with microwave reactors?	73

<b>6.</b>	<b>Applications Examples</b>	<b>74</b>
6.1.	Common Applications	74
6.1.1.	<i>Metal-catalyzed Reactions</i>	74
6.1.2.	<i>Condensation Reactions</i>	77
6.1.3.	<i>Addition Reactions</i>	79
6.1.4.	<i>Substitution Reactions</i>	81
6.1.5.	<i>Rearrangements</i>	83
6.1.6.	<i>Synthesis of and with Ionic Liquids</i>	84
6.1.7.	<i>Nanomaterials</i>	86
6.1.8.	<i>Polymers</i>	90
6.2.	Special Applications	93
6.2.1.	<i>Gaseous Reagents</i>	93
6.2.2.	<i>Chemistry in Water, Near-critical Water, and Supercritical Fluids</i>	94
6.2.3.	<i>Microwave-mediated Synthesis Utilizing Silicon Carbide Tools</i>	97
6.3.	Green Chemistry	103
6.3.1.	<i>General Considerations</i>	103
6.3.2.	<i>Savings under sealed-vessel conditions</i>	105
6.4.	Conclusion	106

<b>7.</b>	<b>Further Reading</b>	<b>108</b>
7.1.	Scientific Publications Involving Anton Paar's Microwave Synthesis Instrumentation	108
7.2.	Recommended Text Books on Microwave Synthesis	124