Anton Paar

SmartRef

Safety Guide

EN	Safety Instructions (Original)	4
DE	Sicherheitshinweise (Übersetzung)	9
FR	Consignes de sécurité (traduction)	14
IT	Istruzioni di sicurezza (traduzione)	19
ES	Instrucciones de seguridad (Traducción)	24
PT	Instruções de segurança (tradução)	29
SV	Säkerhetsanvisningar (översättning)	34
FI	Turvallisuusohjeet (käännös)	39
DA	Sikkerhedsinstruktioner (Oversættelse)	44
NL	Veiligheidsinstructies (vertaling)	49
PL	Instrukcje dotyczące bezpieczeństwa (tłumaczenie)	54
CS	Bezpečnostní pokyny (překlad)	59
SK	Bezpečnostné pokyny (preklad)	64
HU	Biztonsági utasítások (fordítás)	69
HR	Sigurnosne upute (prijevod)	74
SL	Varnostna navodila (prevod)	79
BG	Инструкции за безопасност (Превод)	84
RO	Instrucțiuni de siguranță (Traducere)	89
LT	Saugos nurodymai (vertimas)	94
ET	Ohutusjuhend (tõlge)	99
LV	Drošības norādes (tulkojums)	104
EL	Οδηγίες ασφαλείας (Μετάφραση)	109
ZH	安全说明 (译文)	114

SAFETY INSTRUCTIONS (ORIGINAL)

Anton Paar hereby declares that the radio equipment type SmartRef meets the Directive 2014/53/EU. The full text of the EU Declaration of Conformity and the UK Declaration of Conformity is available at the following Internet address: <https://www.anton-paar.com/easydens#product-documents>

Read this instruction manual before using the device.

Make this instruction manual easily accessible to all persons working with the device.

Follow all tips and instructions in this instruction manual to ensure the correct use and safe functioning of the device.

CONVENTIONS FOR SAFETY MESSAGES

WARNING

- > Warning indicates a hazardous situation which, if not avoided, could result in death or serious injury.

CAUTION

- > Caution indicates a hazardous situation which, if not avoided, could result in minor or moderate injury.

LIABILITY

- > This instruction manual does not claim to address all safety issues associated with the use of the device and samples. It is your responsibility to establish health and safety practices and to determine the applicability of regulatory limitations.
- > Anton Paar warrants the proper functioning of the device only if no modifications are made to mechanics, electronics, or device software.
- > Use the device only for the purpose described in the instruction manual. Anton Paar is not liable for damages caused by incorrect use of the device.
- > The results delivered by the device depend not only on the correct functioning of the device, but also on various other factors.
- > Do not base any decisions concerning safety or commercial issues on data measured by a SmartRef.

INSTALLATION AND USE

WARNING

- > Do not use the device in medical areas.
- > Use the device as a stand-alone unit (the SmartRef cannot be operated without a smartphone or tablet). Do not combine the device with other instruments or devices for automation purposes!
- > Use only accessories and consumables supplied or approved by Anton Paar.
- > Ensure that all operators are familiar with the safety instructions for the device.
- > Keep the device away from children.
- > Ensure that the device is sufficiently supervised during operation.

WARNING

- > In case of damage or malfunction, do not continue operating the device. Do not operate the device under conditions which could result in damage to goods or injuries or loss of life.

OPERATION IN AREAS WITH RISK OF EXPLOSION

WARNING

- > The device is not explosion-proof and therefore must not be operated in areas with risk of explosion.

PERSONAL SAFETY

WARNING

- > Always keep a distance of at least 15 cm (6 in) between the device and an implanted cardiac pacemaker or defibrillator.

CAUTION

- > During operation of the device, the operator must be equipped with appropriate personal protective equipment (safety goggles, gloves).

GENERAL PRECAUTIONS

- > Observe and adhere to your national safety regulations regarding the handling of all substances associated with your measurements (e.g. use safety goggles, gloves, respiratory protection, etc.).
- > For the handling of substances, abide by the corresponding material safety data sheets.

WARNING

- > Fill only substances which will not pose a risk if they leak from the device in case of damage or malfunction.

WARNING

- > Before a measurement, check the wetted parts of the device for chemical resistance to the samples and cleaning agents used.
- > Fill only liquids which will not attack the wetted parts of the device.

WARNING

- > Take care that the liquids (samples and cleaning agents) that you use are chemically compatible when they come into contact with each other. They must not react exothermally or produce harmful substances.
- > Take measures that spilled liquids cannot get into plug connections or venting slots of electrical appliances.
- > Do not leave the device in a location exposed to direct sunlight or near a heat source for any extended period of time.
- > Do not change the ambient temperature of the device abruptly.

CAUTION

- > Be careful with hot samples. Handling samples with temperatures of more than 50 °C (122 °F) bears the risk of burns.

PRECAUTIONS FOR FLAMMABLE SAMPLES AND CLEANING AGENTS

CAUTION

- > Keep potential sources of ignition, like sparks or open flames, at a safe distance from the device.
- > Store only the minimum required amount of sample, cleaning agents, and other flammable materials near the device.
- > Do not spill sample/cleaning agents or leave their containers uncovered. Immediately remove spilled sample/cleaning agents.
- > Ensure that the setup location is sufficiently ventilated. The environment of the device must be kept free from flammable gases and vapors.
- > Provide fire-extinguishing equipment.

BATTERY HANDLING

- > Make sure you insert the batteries with the correct polarity.
- > Keep batteries away from children.
- > Do not leave batteries lying around unattended. They can be dangerous for children or pets if they use them as toys.
- > If a battery has been swallowed, seek medical advice immediately.
- > Leaking or damaged batteries can cause burns if they come into contact with skin. Use gloves for their handling.
- > Never short-circuit or open batteries.
- > Do not expose batteries to heat or throw them into fire.
- > Do not charge non-rechargeable batteries. There is a risk of explosion.

WARNING

- > Do not insert damaged batteries in the battery compartment. They can cause permanent damage to the device and fire.

WARNING

- > Do not use lithium ion rechargeable batteries or lithium ion batteries with the device.
- > Use only batteries type AAA LR03 (alkaline) or rechargeable batteries type AAA HR03 (NiMH).

**WARNING**

- > In case you are not using the device for a long time, remove the batteries. Batteries could leak or be damaged.

**CAUTION**

- > Do not submerge the device in water or any cleaning agents or other liquids.

**CAUTION**

- > Turn the device off prior to exchanging the batteries.
- > Make sure the sample area is empty and dry before exchanging the batteries to avoid spillage of liquids.
- > Only open the device in a clean and dry environment.
- > The inside of the device does not require any cleaning. Do not use liquid cleaning agents on interior parts of the device.
- > Do not open the compartment containing the optical and electrical components of the device.

DEVICE RETURNS

- > For returns, contact the vendor of the device. The device must be cleaned before return.
- > You must not return devices which are contaminated by radioactive materials, infectious agents, or other harmful substances that can pose health hazards.

DISPOSAL

- > Concerning the disposal of the device, observe the legal requirements in your country.

Address of the device producer:
Anton Paar ConsumerTec GmbH
 Kärntner Straße 418
 A-8054 Graz / Austria – Europe

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