**Anton Paar**

# **EasyDens**

## Safety Guide

EN	Safety Instructions (Original)	4
DE	Sicherheitshinweise (Übersetzung)	9
FR	Consignes de sécurité (traduction)	14
IT	Istruzioni di sicurezza (traduzione)	19
ES	Instrucciones de seguridad (Traducción)	24
PT	Instruções de segurança (tradução)	29
SV	Säkerhetsanvisningar (översättning)	34
FI	Turvallisuusohjeet (käännös)	39
DA	Sikkerhedsinstruktioner (Oversættelse)	44
NL	Veiligheidsinstructies (vertaling)	49
PL	Instrukcje dotyczące bezpieczeństwa (tłumaczenie)	54
CS	Bezpečnostní pokyny (překlad)	59
SK	Bezpečnostné pokyny (preklad)	64
HU	Biztonsági utasítások (fordítás)	69
HR	Sigurnosne upute (prijevod)	74
SL	Varnostna navodila (prevod)	79
BG	Инструкции за безопасност (Превод)	84
RO	Instrucțiuni de siguranță (Traducere)	89
LT	Saugos nurodymai (vertimas)	94
ET	Ohutusjuhend (tõlge)	99
LV	Drošības norādes (tulkojums)	104
EL	Οδηγίες ασφαλείας (Μετάφραση)	109
ZH	安全说明 (译文)	114

## SAFETY INSTRUCTIONS (ORIGINAL)

Anton Paar GmbH hereby declares that the radio equipment type EasyDens meets the Directive 2014/53/EU. The full text of the EU Declaration of Conformity and the UK Declaration of Conformity is available at the following Internet address: <https://www.anton-paar.com/corp-en/services-support/document-finder/>

Read this instruction manual before using the instrument.

Make this instruction manual easily accessible to all persons working with the instrument.

Follow all tips and instructions in this instruction manual to ensure the correct use and safe functioning of the instrument.

Warning indicates a hazardous situation which, if not avoided, could result in death or serious injury.

### LIABILITY

- > This instruction manual does not claim to address all safety issues associated with the use of the instrument and samples. It is your responsibility to establish health and safety practices and to determine the applicability of regulatory limitations.
- > Anton Paar GmbH warrants the proper functioning of the instrument only if no modifications are made to mechanics, electronics, or instrument software.
- > Use the instrument only for the purpose described in the instruction manual. Anton Paar GmbH is not liable for damages caused by incorrect use of the instrument.
- > The results delivered by the instrument depend not only on the correct functioning of the instrument, but also on various other factors.
- > Do not base any decisions concerning safety or commercial issues on data measured by EasyDens.

### INSTALLATION AND USE

#### WARNING!

- > Do not use the instrument in medical areas.
- > Use EasyDens only as a stand-alone unit. (The EasyDens cannot be operated without a smart phone or tablet. Therefore, the smartphone is considered as a part of the measurement unit). Do not combine the instrument with other instruments or devices for automation purposes.
- > The instrument is intended for indoor use only.
- > Use only accessories and consumables supplied or approved by Anton Paar GmbH.
- > Ensure that all operators are familiar with the safety instructions for the instrument.
- > Keep the device away from children.
- > Ensure that the instrument is sufficiently supervised during operation.

#### WARNING!

- > In case of damage or malfunction, do not continue operating the instrument. Do not operate the instrument under conditions which could result in damage to goods or injuries or loss of life.
- > Do not expose the instrument to temperatures below 0 °C (32 °F) if the measuring cell contains water. Freezing water will cause rupture of the measuring cell.

#### WARNING!

- > Do not apply any overpressure to the measuring cell. Fill the measuring cell by hand using syringes only
- > Liquids may spurt out of the instrument. Always use the instrument with a waste hose.

#### WARNING! RISK OF EYE INJURIES

- > Liquids may spurt out of the instrument. Always use the instrument with a waste hose. Also observe the precautions found in the Personal Safety section and wear safety goggles and protective gloves

### OPERATION IN AREAS WITH RISK OF EXPLOSION

#### WARNING!

- > The instrument is not explosion-proof and therefore must not be operated in areas with risk of explosion.

### PERSONAL SAFETY

- > Always keep a distance of at least 15 cm (6 in) between the instrument and an implanted cardiac pacemaker or defibrillator.
- > During operation of the instrument, the operator must be equipped with appropriate protective clothing (safety goggles, gloves).

### GENERAL PRECAUTIONS

- > Observe and adhere to your national safety regulations regarding the handling of all substances associated with your measurements (e.g. use safety goggles, gloves, respiratory protection, etc.).
- > For the handling of substances, abide by the corresponding material safety data sheets.

**WARNING!**

- Fill only substances which will not pose a risk if they leak from the instrument in case of damage or malfunction.

**WARNING!**

- Before a measurement, check the wetted parts of the instrument for chemical resistance to the samples and cleaning agents used.
- Fill only liquids which will not attack the wetted parts of the instrument (see section wetted parts).

**WARNING!**

- Take care that the liquids (samples and cleaning agents) that you use are chemically compatible when they come into contact with each other. They must not react exothermally or produce harmful substances.
- Before you start a measurement or cleaning procedure, take care that all parts, in particular the measuring cell, the syringe and the hose for waste are properly connected and in good condition.
- Before you start a measurement or cleaning procedure, check the connection between syringe and the hose for leak tightness.
- Take measures that spilled liquids cannot get into plug connections or venting slots of electrical appliances.
- Do not leave the instrument in a location exposed to direct sunlight or near a heat source for any extended period of time.
- Do not change the ambient temperature of the instrument suddenly.

**WARNING!**

- Do not expose the instrument to temperatures below 0 °C (32 °F) if the measuring cell contains water. Freezing water will cause rupture of the measuring cell.

**HOT SAMPLES**

- Be careful with hot samples. Handling samples with temperatures of more than 50 °C (122 °F) bears the risk of burns.
- Observe the maximum sample temperature. At temperatures over 80 °C (176 °F), plastic parts soften and connections become leaky.

**PRECAUTIONS FOR FLAMMABLE SAMPLES AND CLEANING AGENTS**

- Keep potential sources of ignition, like sparks or open flames, at a safe distance from the instrument.
- Store only the minimum required amount of sample, cleaning agents, and other flammable materials near the instrument.
- Do not spill sample/cleaning agents or leave their containers uncovered. Immediately remove spilled sample/cleaning agents.

**WARNING!**

- If the U-tube breaks, wipe off leaked liquids immediately to avoid risk of fire.
- Ensure that the setup location is sufficiently ventilated. The environment of the instrument must be kept free of flammable gases and vapors.
- Provide fire-extinguishing equipment.

**BATTERY HANDLING**

- Please see how to insert the batteries into the device as displayed in picture - View of how to add batteries and correct polarity. Make sure you put the batteries in according to the correct polarity.
- Keep batteries away from children.
- Do not leave batteries lying around unattended. They can be dangerous for children or pets if they use them as toys.
- If a battery has been swallowed, seek medical advice immediately.
- Leaking or damaged batteries can cause burns if they come into contact with your skin. Use gloves for their handling.
- Never short-circuit or open batteries.
- Do not expose batteries to heat or throw them into fire.
- Do not charge non-rechargeable batteries. There is a risk of explosion.

**WARNING!**

- Do not insert damaged batteries in the battery compartment. They can cause short circuits and fire.

**WARNING!**

- Do not use lithium ion rechargeable batteries or lithium ion batteries with the instrument.
- Use only batteries type AA LR6 (alkaline) or rechargeable batteries type AA HR6 (NiMH).

- > When replacing batteries, this step should be performed in an environment where there is relatively low humidity

### WARNING!

- > In the case that you are not using the instrument for a long time, remove the batteries. Batteries could leak or be damaged.

### INSTRUMENT RETURNS

- > For returns, contact the vendor of the instrument. The instrument must be cleaned before return.
- > You must not return instruments which are contaminated by radioactive materials, infectious agents, or other harmful substances that cause health hazards.

### DISPOSAL

- > Concerning the disposal of the instrument, observe the legal requirements in your country.

Address of the instrument producer:  
**Anton Paar** GmbH  
 Anton-Paar-Str. 20  
 A-8054 Graz/Austria – Europa

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